

Amsco Autoclave Operation and Programming Instructions

A. To run:

1. Look at the screen. It will show you which programs are for a liquid cycle (slow exhaust), and which are gravity (solids, fast exhaust).
2. Press the number of the program that you desire (#1-4). The screen will then show additional parameters of the program (times, etc.). You may have to browse through several until you find one that you like. Don't assume that a program that you used last time is still the same!! If you need to make a new program, see Section B.
3. To start, press the number of the program a second time. if you wait too long, the display will reset, and you will have to go back to step 2.

B. To change program parameters:

1. Lift up the sliding panel.
2. Press "change values". Wait until the date and time screen appear before doing anything else.
3. Press "change values" again. This will cause the screen for program #1 to appear. If you want to change program #1, skip to step 4. If you want to change another program, keep pressing "change values" until that program appears.
4. Press the horizontal arrows to scroll to the parameter that you want to change.
5. Press the vertical arrows to change the values (such as time).
6. When done making the changes, press "save values."
7. Once the printer has completed printing the new program parameter list (you have to wait), then start the program as described in Section A2-3 (in other words, just press the number of the program twice).